



NatuTec GmbH Forschungs- und Laborsysteme
 Königslacher Str. 15-21
 60528 Frankfurt am Main
 phone: +49 (0)69 677 378-0
 fax: +49 (0)69 677 378-79
 email: info@natutec.com
 http://www.natutec.com

Taq DNA Polymerase

Cat. No.:	TP-0200-A1	200 u Taq DNA Polymerase (5 u/μl), buffer A with MgCl₂
	TP-1000-A1	1000 u Taq DNA Polymerase (5 u/μl), buffer A with MgCl₂
	TP-0200-A0	200 u Taq DNA Polymerase (5 u/μl), buffer A without MgCl₂
	TP-1000-A0	1000 u Taq DNA Polymerase (5 u/μl), buffer A without MgCl₂
	TP-0200-B1	200 u Taq DNA Polymerase (5 u/μl), buffer B with MgCl₂
	TP-1000-B1	1000 u Taq DNA Polymerase (5 u/μl), buffer B with MgCl₂
	TP-0200-B0	200 u Taq DNA Polymerase (5 u/μl), buffer B without MgCl₂
	TP-1000-B0	1000 u Taq DNA Polymerase (5 u/μl), buffer B without MgCl₂

Source: Isolated from a recombinant *E. coli* strain.

Description: Taq DNA polymerase is a thermostable enzyme of approximately 94 kDa isolated from eubacterium *Thermus aquaticus* strain YT-1. This unmodified enzyme replicates DNA at 72 °C. The enzyme catalyzes the polymerization of nucleotides into duplex DNA in the 5'-3' direction in the presence of magnesium ions and possesses a 5'-3' exonuclease activity. The enzyme is highly purified and is free of nonspecific endo- or exonucleases. Taq DNA polymerase adds extra 3'-dA nucleotide(s) overhangs to its reaction products.

Concentration: 5 u/μl

Storage conditions: 20 mM Tris-HCl (pH 7.6); 100 mM KCl; 0.1 mM EDTA; 1 mM DTT; 0.5% Triton X-100; 50% glycerol.
Store at -20 °C.

Reaction buffer A w MgCl₂ (1x): 60 mM Tris-HCl (pH 8.5 at 25 °C); 1.5 mM MgCl₂; 25 mM KCl; 10 mM 2-mercaptoethanol; 0.1% Triton X-100

Reaction buffer A w/o MgCl₂ (1x): 60 mM Tris-HCl (pH 8.5 at 25 °C); 25 mM KCl; 10 mM 2-mercaptoethanol; 0.1% Triton X-100

Reaction buffer B w MgCl₂ (1x): 16 mM (NH₄)₂SO₄; 67 mM Tris-HCl (pH 8.8); 0.1% Tween 20; 2.5 mM MgCl₂

Reaction buffer B w/o MgCl₂ (1x): 16 mM (NH₄)₂SO₄; 67 mM Tris-HCl (pH 8.8); 0.1% Tween 20

Unit definition: One unit of enzyme catalyzes the incorporation of 10 nanomoles of dNTPs into adsorbable form in 30 minutes at 72 °C.

Units assay conditions: 60 mM Tris-HCl (pH 8.5); 25 mM KCl; 1.5 mM MgCl₂; 0.1% Triton X-100; 10 mM 2-mercaptoethanol; 200 μM dATP, dCTP, dGTP; 50 μM ³H-TTP; 12.5 μg activated calf thymus DNA.

Reaction Conditions (for a 50 µl volume)

10x Taq DNA Polymerase buffer	5 µl
MgCl ₂	as required
40 mM dNTP Mix	0.5 – 1.0 µl
Primer I (0.1 - 1 µM)	as required
Primer II (0.1 - 1 µM)	as required
Template DNA (10 pg - 1 µg)	as required
NatuTec Taq DNA Polymerase (5 u/µl)	0.25 – 1.0 µl
(50 mM MgCl ₂ Solution)	(1.5 – 4.0 µl)
Water (ddH ₂ O)	up to 50 µl


Initial Denaturing: 94 – 96°C for 5 min
Denaturing: 94 – 96°C for 45 sec
Annealing: depends on your template
Elongation : 70 – 72°C (allowing 30 – 60 sec/kb)
Final Elongation: 70 – 72°C for 7 min
Cycles: 25 – 30

6 kb amplification control:

Template was genomic DNA of *Staphylococcus aureus*.

Lane 1: 1 kb DNA marker

Lane 2: amplified 6 kb DNA fragment

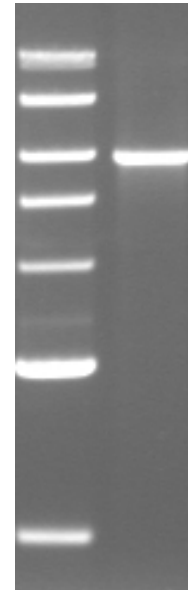

Please note:

This data is intended for use as a guide only.
Conditions will vary from reaction to reaction and may need optimisation.

Patent Disclaimer:

Unless explicitly stated, no licence or immunity under any patent is either granted or implied by the sale of any of our products. *NatuTec GmbH* does not warrant that the resale or use of its products delivered will not infringe the claims of any patents, trademarks or copyright covering use of the product itself, or its use in combination with other any other products, or its use in the operation of any process. Furthermore, the purchaser assumes all risks of patent, trademark or copyright infringement associated with any such use, combination or operation.

Einige Anwendungen, für die dieses Produkt eingesetzt werden kann, sind in bestimmten Ländern patentrechtlich geschützt. Da mit seinem Kauf keine Lizenzen für die Verwendung im Rahmen patentrechtlich geschützter Anwendungen erworben werden, kann, abhängig von der Verwendung und abhängig vom Anwendungsland, der Erwerb entsprechender Lizenzen erforderlich sein.